



BRT test tubes – instruction manual heating block

Start of operation

- Connect power cord to 230 VAC mains
- Turn on „Netz“ main switch, yellow LED „Heizen“ is on
- After 10 minutes the yellow LED „Heizen“ starts flashing and indicates that the heating bloc is operable and temperature has reached 65.0 °C

Specification

Temperature: 65.0 °C +/- 0.3 °C
Voltage: 230 VAC, 30 watt

Temperature check

- Use appropriate thermometer, e.g. digital thermometer AiM GmbH art. no. 3057
- Put the thermometer in the small drilling and read temperature after a few minutes.

Attention

- Safeguard heating block against crushing, water and moisture.
- Dry cleaning.
- We recommend operating the heating bloc for inhibitor testing at room temperature.

